

MINLITE SERIES

When your **Minilite** laser arrives, contact Continuum immediately if there is damage to the crates or boxes. If no damage is evident, unpack and move the laser system to your lab, consulting the Installation section of your laser manual for details. After unpacking, again check for visible damage. The **Minilite** lasers are customer installed. However, if you would like to receive training and have your laser installed by one of our technicians, please call our Service Manager at 408-727-3240 (ask for service) to receive a quote.

We suggest that you have the equipment and materials listed below available during installation. These items will also be useful for operating and maintaining your system later.

- 200-220V single phase power, 8 amp service, and a plug to fit your outlet
- Laser safety glasses
- Methanol, 50ml, reagent grade or better
- Distilled water, 1.5 gallons (6 liters)
- Kentek zap It burn paper
- Power meter, 10 Watt, with volume absorbing detector
- Mini Mag Lite and Inspection mirror (dental type)
- IR phosphor card
- Lens tissue paper

If you have any questions or would like additional information, please call our Technical Support Department at (877)272-7783.
