

SURELITE III

When your Surelite III laser system arrives, contact Continuum immediately if there is damage to the crates or boxes. If no damage is evident, unpack and move the laser system to your lab, consulting the Installation section of your laser manual for details. After unpacking, again check for visible damage and call Continuum to schedule your installation.

Continuum asks that you have the equipment and materials listed below available during installation. These items will also be useful for operating and maintaining your system later. Please contact us in the Service Department of our Santa Clara office if you cannot provide all of them.

- Laser safety glasses
- Methanol, 50 ml, reagents grades or better
- Distilled water, 1.5 gallons (6 liters)
- 200-220 V single phase power, 8 amp service, and a plug to fit your outlet
- Kentek Zap It Burn Paper
- Power meter, 10 watt, with volume absorbing detector
- Mini Mag Lite and Inspection mirror (dental type)
- IR phosphor card
- Lens tissue paper

If you have any questions or would like additional information, please don't hesitate to call our Technical Support Dept @ (877)272-7783.
